


WellMate Setup Instructions for 6, 12, 24, & 48 well Plate Dispensing

6-Well Plate Set-Up

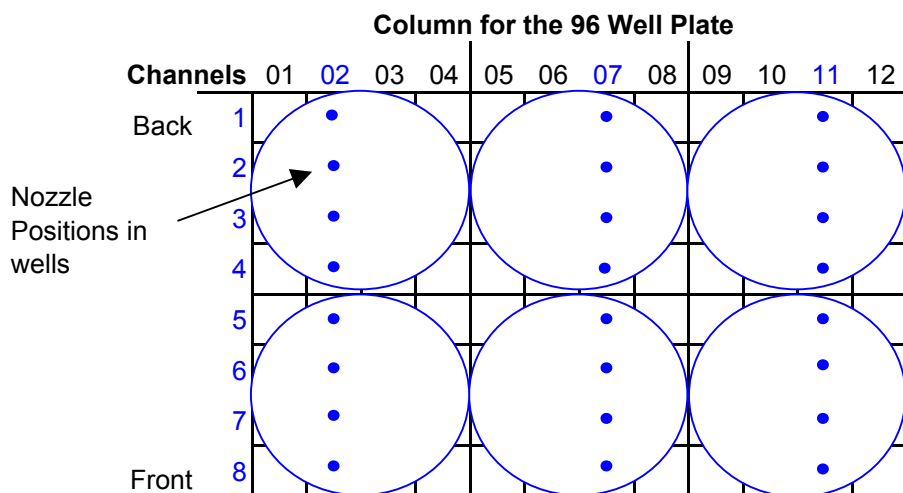


- Indicates location where fluid is dispensed in well
Set volume @ 1/4 of desired amount
For Example: Set volume @500uL if desired amount is 2000uL
Leave nozzle arm in 96 channel mode

How to program the WellMate for 6 Well Plates

Activated Columns	
Column Number	ON(1)/OFF(0)
01	0
02	1
03	0
04	0
05	0
06	0
07	1
08	0
09	0
10	0
11	1
12	0

Activated Channels	
Channels	Connected/Disconnected
Back	1 C
	2 C
	3 C
	4 C
	5 C
	6 C
	7 C
Front	8 C

12-Well Plate Set-Up

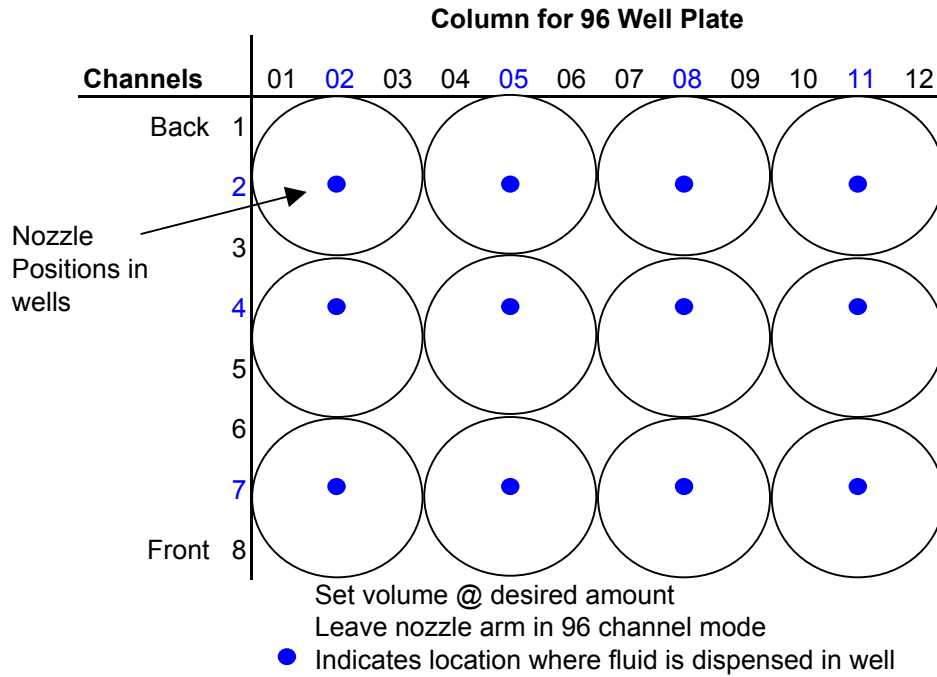


How to program the WellMate for 12 Well Plates

Activated Columns	
Column Number	ON(1)/OFF(0)
01	0
02	1
03	0
04	0
05	1
06	0
07	0
08	1
09	0
10	0
11	1
12	0

Activated Channels	
Channels	Connected/Disconnected
Back 1	D
2	C
3	D
4	C
5	D
6	D
7	C
Front 8	D

24-Well Plate Set-Up

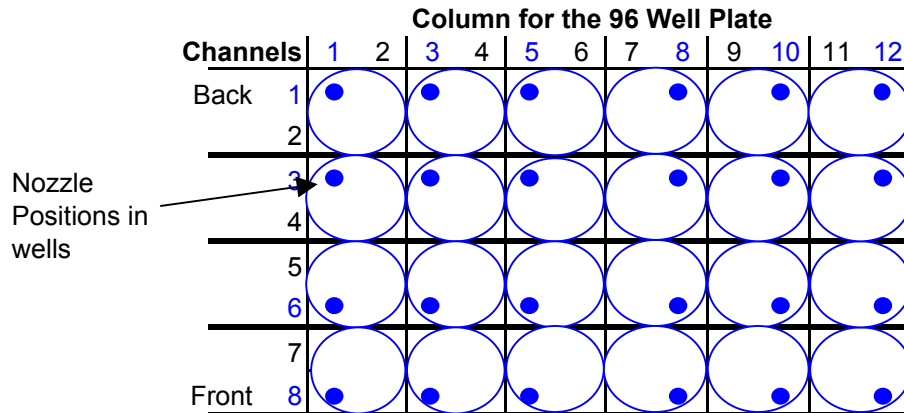

Set volume @ desired amount

Leave nozzle arm in 96 channel mode


- Indicates location where fluid is dispensed in well

How to program the WellMate for 24 Well Plates

Activated Columns	
Column Number	ON(1)/OFF(0)
01	1
02	0
03	1
04	0
05	1
06	0
07	0
08	1
09	0
10	1
11	0
12	1

Activated Channels	
Channels	Connected/Disconnected
Back 1	C
2	D
3	C
4	D
5	D
6	C
7	D
Front 8	C

48-Well Plate Set-Up

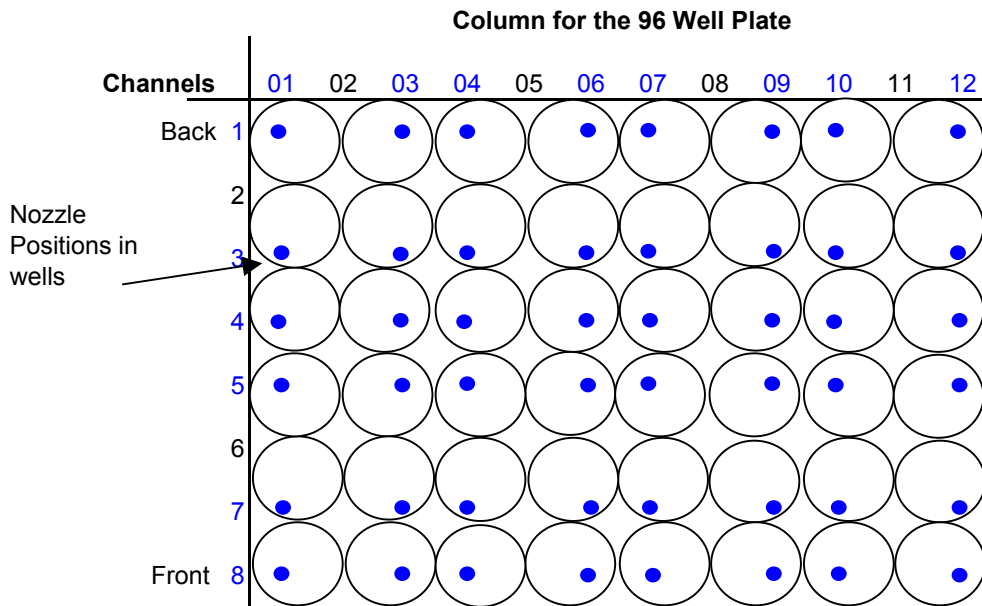

- Set volume @ desired amount
- Leave nozzle arm in 96 channel mode
- Indicates location where fluid is dispensed in well

How to program the WellMate for 48 Well Plates

Activated Columns	
Column Number	ON(1)/OFF(0)
01	1
02	0
03	1
04	1
05	0
06	1
07	1
08	0
09	1
10	1
11	0
12	1

Activated Channels	
Channels	Connected/Disconnected
Back	1 C
	2 D
	3 C
	4 C
	5 C
	6 D
	7 C
Front	8 C